




LABORATORY SECTOR

AUTOCLAVE



Kalstein®

A different accompaniment, at your service

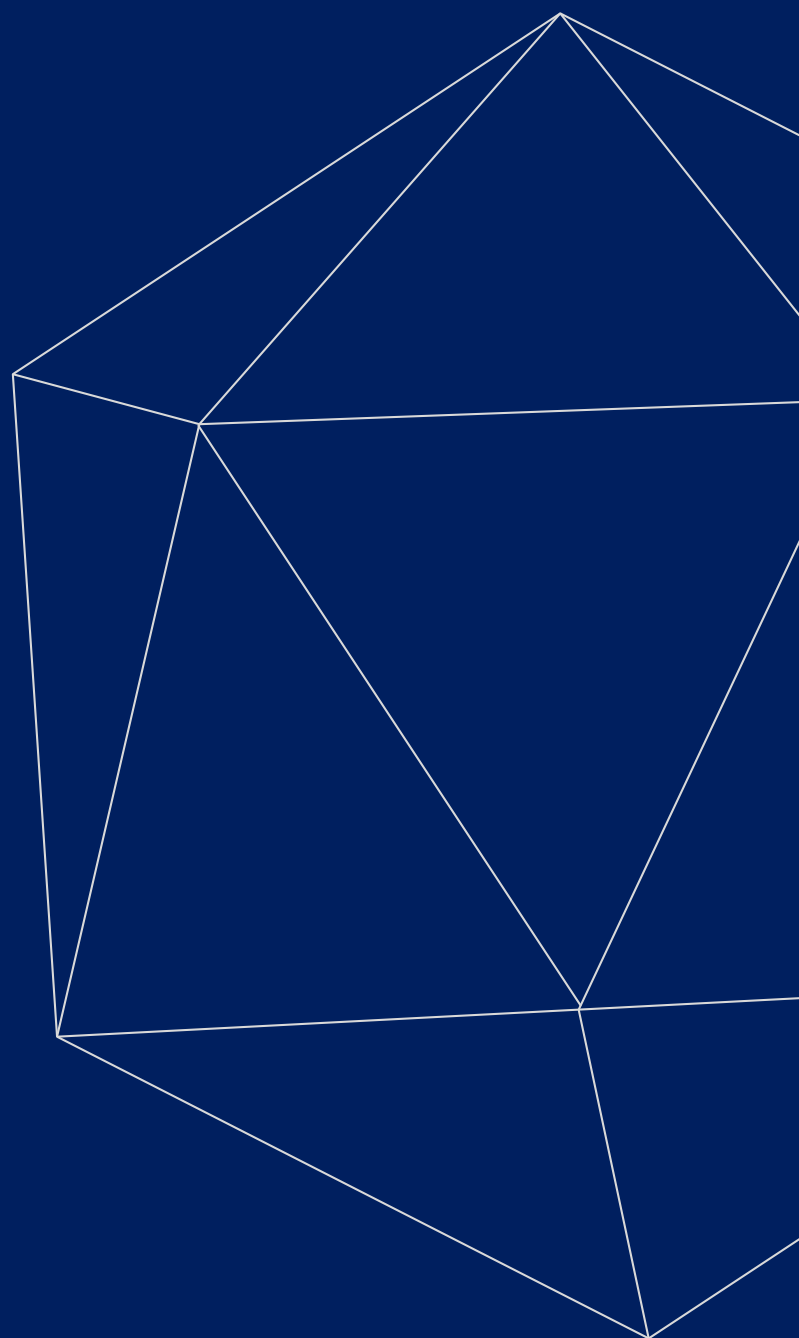



Kalstein®

A different accompaniment, at your service

OUR SERVICES

Benefits and Support

In Kalstein France, we take care of the full satisfaction of our customers, that is why we provide value-added services of the highest level based on our experience.



Online Inductions and Trainings

In any part of the world, receive your induction or training from our specialized team of engineers



Quick Response

Our work team is always available to response all your consults or questions, in order to support you in any situation.



#Letsgivemore

Thanks to your purchase, a donation will be made to a non-profit foundation that fights against breast cancer and helps most vulnerable communities.



Technical Support

Enjoy of personalized advice for the correct preventive and corrective maintenance of your equipment, thanks to Kalstein's manuals and articles, special catalogues and video tutorials.



Delivery Logistics

We take care of all the necessary logistics for the dispatch of your goods, whether is by sea, land or air.



Kalstein Worldwide

With more than 25 years growing with our customers, Kalstein's multiformat and modern content, is now present in more than 10 countries and increasing.



KALSTEIN® Autoclave
Class B Benchtop Autoclave
Model YR03385B // YR03385C



#Letsgivemore ♥

Thanks to your purchase, a donation will be made to a non-profit foundation that fights against breast cancer and helps most vulnerable communities.



All rights reserved © KALSTEIN France S. A. S.,
2 Rue Jean Lantier • 75001 Paris •
+33 1 78 95 87 89 / +33 6 80 76 07 10 •
<https://kalstein.eu>
KALSTEIN FRANCE, S. A. S



KALSTEIN® Autoclave Class B Benchtop Autoclave Model YR03385B // YR03385C



Product Description:

8 liter, 12 liter, 18 liter and 23 liter tabletop autoclaves with class B pre-vacuum, complies with the European standard EN13060. These autoclaves require minimal space, being an ideal solution for small businesses such as tattoo, podiatric consultations, beauty salons, veterinary medicine and laboratories.

Excellent Performance:

- ✓ Fully automatic tabletop autoclaves with pre-programmed sterilization cycles.
- ✓ Chamber and door in stainless steel, with electro-polished finish.
- ✓ High precision control system for perfect sterilization results.
- ✓ Independent steam generator, provides enough steam for faster cycles.
- ✓ Friendly interface for the user, it can be easily operated by the user.
- ✓ LCD screen, continuously displays information about the cycle process.
- ✓ Open type water tank, this feature makes it easier to fill water and clean the tank.

Monitoring and Security:

- ✓ The cycle will not start unless the door is properly closed.
- ✓ Automated system with double safety locking device, prevents the door from opening while the chamber is pressurized (optional).
- ✓ Safety valves prevent generator and chamber overpressure.
- ✓ A safety thermostat prevents overheating in the chamber and the generator.
- ✓ An automatic safety lock prevents accidental opening of the door during the sterilization process.
- ✓ An electricity protector automatically cuts the power if somewhere you have a short circuit or the power supplier is not very stable.
- ✓ Failure alarm, indicates a cycle failure and details the error code.
- ✓ Door alarm, indicates that the door is not secured.
- ✓ Low water alarm, indicates when the water tank is empty or too full.

Documentation:

- ✓ Internal Printer: prints the detailed information of the program cycle. (optional).
- ✓ USB port and USB memory: all the information of the program cycle is automatically described on the USB memory and can be read directly on any PC. (optional)
- ✓ Internal memory: the information of the last 20 cycles is automatically stored in the autoclave system, which can be printed at any time.



#Letsgivemore

Thanks to your purchase, a donation will be made to a non-profit foundation that fights against breast cancer and helps most vulnerable communities.



All rights reserved © KALSTEIN France S. A. S.,
2 Rue Jean Lantier • 75001 Paris •
+33 1 78 95 87 89 / +33 6 80 76 07 10 •
<https://kalstein.eu>
KALSTEIN FRANCE, S. A. S



KALSTEIN® Autoclave
Class B Benchtop Autoclave
Model YR03385B // YR03385C



Technical Description:

Order Code	YR03385B	YR03385C
Chamber Dimension	Ø319×417 mm	Ø319×617 mm
Chamber Volume	29 L	45 L
Number of Trays	3	3
Material of Trays	Stainless steel	Stainless steel
Size of Trays	240×340 mm	240×540 mm
Voltage, Freq.	220/110V, 50/60Hz	220/110V, 50/60Hz
Power	2300 w	3800 w
Overall Dimension (WxHxD)	640x640x560 mm	840x640x560 mm
Weight	95 kgs	115 kgs



#Letsgivemore ♥

Thanks to your purchase, a donation will be made to a non-profit foundation that fights against breast cancer and helps most vulnerable communities.



All rights reserved © KALSTEIN France S. A. S.,
 2 Rue Jean Lantier • 75001 Paris •
 +33 1 78 95 87 89 / +33 6 80 76 07 10 •
<https://kalstein.eu>
KALSTEIN FRANCE, S. A. S



KALSTEIN® Autoclave
Class B Benchtop Autoclave
Model YR03385B // YR03385C



Technical Description:

Program	Temp. [°C]	Pressure [Mpa]	Sterilization time [min.]
QUICK	134	210	4
WRAPPED	134	210	10
	121	110	30
PRION	134	210	18
DRY	–	–	–
B&D TEST	134	210	3.5
HELIX TEST	134	210	3.5
VACUUM TEST	–	–	–



#Letsgivemore ♥

Thanks to your purchase, a donation will be made to a non-profit foundation that fights against breast cancer and helps most vulnerable communities.



All rights reserved © KALSTEIN France S. A. S.,
 2 Rue Jean Lantier • 75001 Paris •
 +33 1 78 95 87 89 / +33 6 80 76 07 10 •
<https://kalstein.eu>
KALSTEIN FRANCE, S. A. S





 **Kalstein**[®]
A different accompaniment, at your service

#LETSGIVEMORE 

**WITH THE
ACQUISITION
OF A KALSTEIN
EQUIPMENT**

YOU MAKE A CONTRIBUTION TO:



**Fundación
Jacinto Convit**

<https://www.jacintoconvit.org/>



<https://onetreepanted.org>



Humatem

Solidarité et coopération
biomédicale


<https://www.humatem.org/>



**MANIAPURE
FOUNDATION USA**

<https://www.maniapurefoundation.org/>

Articles Kalstein
(Social Networks and Official Websites)

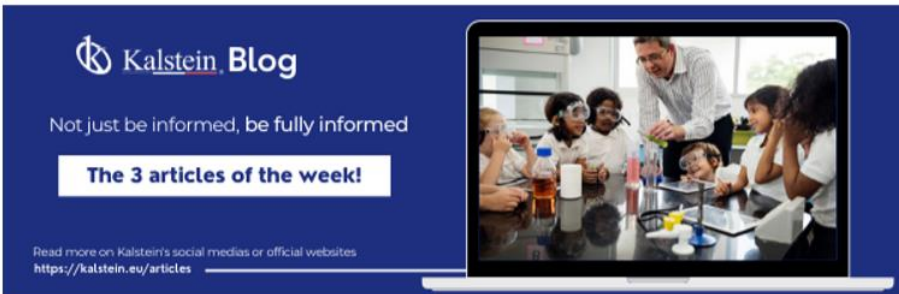

Kalstein Blog
Not just be informed, be fully informed

The 3 articles of the week!

Read more on Kalstein's social medias or official websites
<https://kalstein.eu/articles>



1
Read this ARTICLE NOW

How does the climate affect your microscope?

Kalstein Blog

Article of the week:
How does the climate affect your microscope?

Read more on Kalstein's social medias or official websites
<https://kalstein.eu/articles>



2
Read this ARTICLE NOW

What is the importance of the laboratory refrigerator?

Kalstein Blog

Article of the week:
What is the importance of the laboratory refrigerator?

Read more on Kalstein's social medias or official websites
<https://kalstein.eu/articles>

Kalstein Training Videos
(Youtube Channel and Official Websites)



Kalstein France - Jacinto Convit Foundation / Allianc...

23 views • 3 years ago



Kalstein's Microscopes

445 views • 1 week ago



PARIS - FRANCE
HEADQUARTER

2 Rue Jean Lantier
75001 Paris - France

Tlf: +33 (0) 01 78 95 87 89
Cel: +33 (0) 6 80 76 07 10
sales@kalstein.eu

<https://kalstein.eu/>
<https://www.kalstein.fr/>

