



LABORATORY SECTOR

AUTOCLAVE



Kalstein®

A different accompaniment, at your service





Kalstein®

A different accompaniment, at your service

OUR SERVICES

Benefits and Support

In Kalstein France, we take care of the full satisfaction of our customers, that is why we provide value-added services of the highest level based on our experience.



Online Inductions and Trainings

In any part of the world, receive your induction or training from our specialized team of engineers



Quick Response

Our work team is always available to response all your consults or questions, in order to support you in any situation.



#Letsgivemore

Thanks to your purchase, a donation will be made to a non-profit foundation that fights against breast cancer and helps most vulnerable communities.



Technical Support

Enjoy of personalized advice for the correct preventive and corrective maintenance of your equipment, thanks to Kalstein's manuals and articles, special catalogues and video tutorials.



Delivery Logistics

We take care of all the necessary logistics for the dispatch of your goods, whether is by sea, land or air.



Kalstein Worldwide

With more than 25 years growing with our customers, Kalstein's multiformat and modern content, is now present in more than 10 countries and increasing.



KALSTEIN® Autoclave
Tabletop Autoclave Class B
Model YR03387 // YR03390



#Letsgivemore ♥

Thanks to your purchase, a donation will be made to a non-profit foundation that fights against breast cancer and helps most vulnerable communities.



All rights reserved © KALSTEIN France S. A. S.,
2 Rue Jean Lantier • 75001 Paris •
+33 1 78 95 87 89 / +33 6 80 76 07 10 •
<https://kalstein.eu>
KALSTEIN FRANCE, S. A. S



KALSTEIN® Autoclave Tabletop Autoclave Class B Model YR03387 // YR03390



Product Description:

Fully automatic tabletop autoclave, triple fractionated pre-vacuum, works with a precision method controlling pressure and temperature. 8 Liter, 12 Liter, 18 Liter and 23 Liter tabletop autoclaves with class B pre-vacuum, complies with the European standard EN13060. These autoclaves require minimal space, being an ideal solution for small businesses such as tattoo, podiatric consultations, beauty salons, veterinarians and laboratories..

Excellent Performance:

- ✓ Fully automatic tabletop autoclaves with pre-programmed sterilization cycles.
- ✓ Chamber and door in stainless steel, with electro-polished finish.
- ✓ High precision control system for perfect sterilization results.
- ✓ Independent steam generator, provides enough steam for faster cycles.
- ✓ Friendly interface for the user, it can be easily operated by the user.
- ✓ LCD screen, continuously displays information about the cycle process.
- ✓ Open type water tank, this feature makes it easier to fill water and clean the tank.

Monitoring and Security:

- ✓ The cycle will not start unless the door is properly closed.
- ✓ Automated system with double safety locking device, prevents the door from opening while the chamber is pressurized (optional).
- ✓ Safety valves prevent generator and chamber overpressure.
- ✓ A safety thermostat prevents overheating in the chamber and the generator.
- ✓ An automatic safety lock prevents accidental opening of the door during the sterilization process.
- ✓ An electricity protector automatically cuts the power if somewhere you have a short circuit or the power supplier is not very stable.
- ✓ Failure alarm, indicates a cycle failure and details the error code.
- ✓ Door alarm, indicates that the door is not secured.
- ✓ Low water alarm, indicates when the water tank is empty or too full.



#Letsgivemore ♥

Thanks to your purchase, a donation will be made to a non-profit foundation that fights against breast cancer and helps most vulnerable communities.



All rights reserved © KALSTEIN France S. A. S.,
2 Rue Jean Lantier • 75001 Paris •
+33 1 78 95 87 89 / +33 6 80 76 07 10 •
<https://kalstein.eu>
KALSTEIN FRANCE, S. A. S



KALSTEIN® Autoclave Tabletop Autoclave Class B Model YR03387 // YR03390



Technical Description:

Model Number	YR03387	YR03388	YR03389	YR03390
Chamber Dimension	Ø170×320 mm	Ø200×360 mm	Ø247×350 mm	Ø247×450 mm
Chamber Volume	8 L	12 L	18 L	23 L
Number of Trays	2/ Max. 3	2/ Max. 3	3/ Max. 5	3/ Max. 5
Material of Trays	Aluminum	Stainless Steel	Aluminum	Aluminum
Size of Trays.	130x265mm	156x295mm	192x285mm	192x380mm
Voltage Freq.	220/110V	220/110V	220/110V	220/110V
	50/60Hz	50/60Hz	50/60Hz	50/60Hz
POWER.r	2000W	2000W	2000W	2200W
Overall Dimension (WxHxD)	420x370x525 mm	420x370x595 mm	490x455x600 mm	490x455x690 mm
Weight	37 kgs	40 kgs	47 kgs	53 kgs



#Letsgivingore

Thanks to your purchase, a donation will be made to a non-profit foundation that fights against breast cancer and helps most vulnerable communities.



All rights reserved © KALSTEIN France S. A. S.,
2 Rue Jean Lantier • 75001 Paris •
+33 1 78 95 87 89 / +33 6 80 76 07 10 •
<https://kalstein.eu>
KALSTEIN FRANCE, S. A. S



KALSTEIN® Autoclave Tabletop Autoclave Class B Model YR03387 // YR03390



Program	Temp. [-C]	Pressure [Mpa]	Sterilization time [min.]	
SOLID	134	210	4	
	121	110	20	
LIQUID	134	210	10	
	121	110	30	
WRAPPED	134	210	10	
	121	110	30	
TEXTILE	134	210	10	
	121	110	30	
PRION	134	210	18	
DRYING	–	–	–	ene-20
B&D TEST	134	210	3.5	
HELIX TEST	134	210	3.5	
VACUUM TEST	–	–	–	15-20



#Letsgivingore

Thanks to your purchase, a donation will be made to a non-profit foundation that fights against breast cancer and helps most vulnerable communities.



All rights reserved © KALSTEIN France S. A. S.,
2 Rue Jean Lantier • 75001 Paris •
+33 1 78 95 87 89 / +33 6 80 76 07 10 •
<https://kalstein.eu>
KALSTEIN FRANCE, S. A. S





 **Kalstein**[®]
A different accompaniment, at your service

#LETSGIVEMORE 

**WITH THE
ACQUISITION
OF A KALSTEIN
EQUIPMENT**

YOU MAKE A CONTRIBUTION TO:



**Fundación
Jacinto Convit**

<https://www.jacintoconvit.org/>



<https://onetreepanted.org>



Humatem

Solidarité et coopération
biomédicale


<https://www.humatem.org/>



**MANIAPURE
FOUNDATION USA**

<https://www.maniapurefoundation.org/>

Articles Kalstein
(Social Networks and Official Websites)

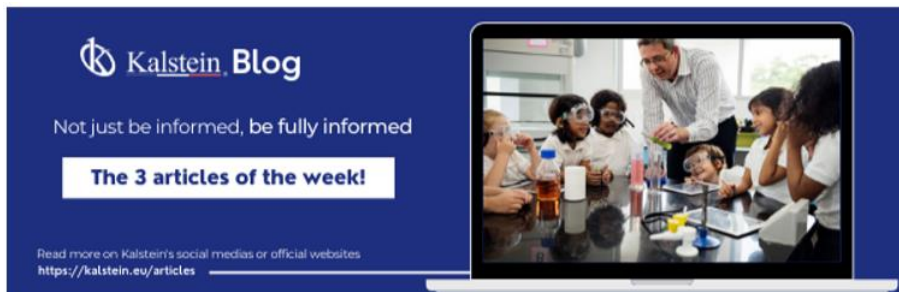

Kalstein Blog
Not just be informed, be fully informed

The 3 articles of the week!

Read more on Kalstein's social medias or official websites
<https://kalstein.eu/articles>



PARIS - FRANCE
HEADQUARTER
2 Rue Jean Lantier
75001 Paris - France

Tlf: +33 (0) 01 78 95 87 89
Cel: +33 (0) 6 80 76 07 10
sales@kalstein.eu

<https://kalstein.eu/>
<https://www.kalstein.fr/>



1
Read this ARTICLE NOW

How does the climate affect your microscope?

Kalstein Blog

Article of the week:
How does the climate affect your microscope?

Read more on Kalstein's social medias or official websites
<https://kalstein.eu/articles>



2
Read this ARTICLE NOW

What is the importance of the laboratory refrigerator?

Kalstein Blog

Article of the week:
What is the importance of the laboratory refrigerator?

Read more on Kalstein's social medias or official websites
<https://kalstein.eu/articles>

Kalstein Training Videos
(Youtube Channel and Official Websites)



Kalstein France - Jacinto Convit Foundation / Allianc...
23 views • 3 years ago



Kalstein's Microscopes
445 views • 1 week ago

